

Attune™ – Daily Maintenance

Beginning of the Day

- 1 POWER ON your system



Notes: if the system is already ON, close and relaunch the software


- 2 Run Startup



- 3 Run Performance Test




- 4 Sanitize Attune SIP

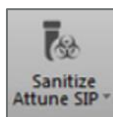


Between Experiments or Plates

- 1 Sanitize SIP with Bleach
- 2 Sanitize SIP with Wash

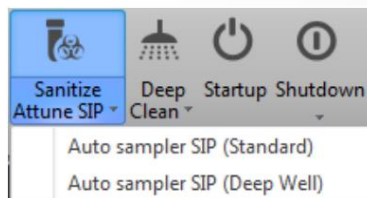


After sticky cells, prefer Deep Clean

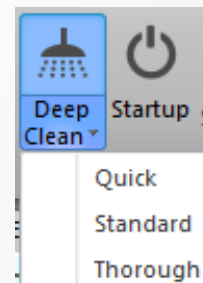

End of the Day

- 1 Sanitize SIP with Flow Cell Cleaning solution diluted 1:3 in Ultra Pure Water (only once a week if <6 hours of use)

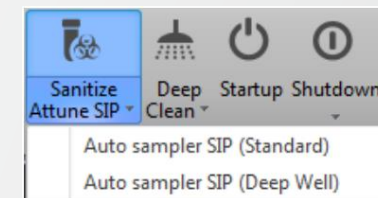

- 2 Run Shutdown

- 3 POWER OFF the entire system at least once a week